

Agar Plates for Phage Inhibition Assay

Add to 1 liter distilled H₂O

| | |
|---------------------------------|-------|
| Bacto-Agar (Stores #436573) | 12 gm |
| Bacto-Tryptone (Stores #436566) | 10 gm |
| NaCl | 8 gm |
| Na Citrate | 2 gm |

Dissolve by heating and stirring.

After above is dissolved, add 3 g glucose.

Auto clave.

Cool till warm to touch and pour ~ 15 ml per plate.

1 liter makes about 70 plates.