

MEDIA FOR PHAGE INHIBITION ASSAY

Soft Agar

Add to 1 liter distilled H₂O

| | |
|----------------|-------|
| Bacto-Agar | 7 gm |
| Bacto-Tryptone | 10 gm |
| NaCl | 8 gm |
| Na Citrate | 2 gm |

Dissolve by heating and stirring. After above is dissolved, add 3 g glucose. Dispense into ~200 ml portions and autoclave.

Nutrient Broth

Add to 1 liter distilled H₂O

| | |
|----------------|------|
| Bacto-Tryptone | 10 g |
| NaCl | 5 g |

Dispense into ~200 ml portions and auto clave.