


Procedure for mouse Ouchterlony

To start fill out an Ouchterlony sheet

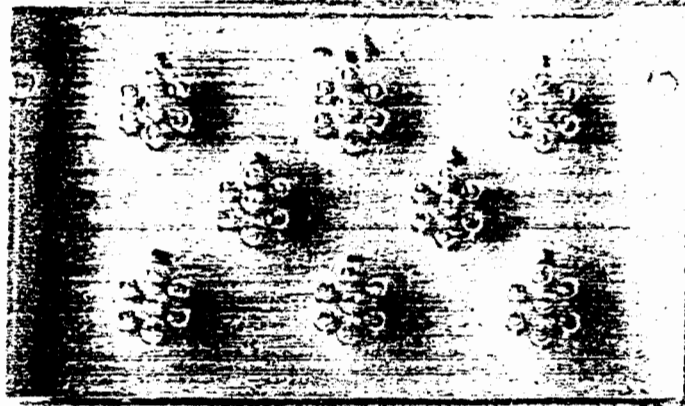

1. Use pre-coated Agar slide
2. Punch slide using the above temp. plate & holder. Suck the agar out with the large needle then clean out wells with the small needle
3. Add serum. Fill wells so that it's eye level with agar.
4. Incubate slide in slide box for 4 hrs. at R.T. ^(or O.N.) _(4°C)
5. wet slide with H₂O, cover w/ bibulous paper and dry for 1 hour.
6. Place slide in 0.5% Sodium Bicarbonate Buffer "Wash" for 4 hours or O.N.
7. Stain slide in Buffalo Black for 3-5 min.
8. Destain in 5% Acetic Acid until background is clear.
9. Dry for 10 min.