

47.

5% Sodium Bicarbonate (10x stock)

For 500 ml:

25 gms. Sodium Bicarbonate  
500 ml d. H<sub>2</sub>O

1. Dissolve sodium Bicarbonate  
in d. H<sub>2</sub>O.

0.5% Sodium Bicarbonate → N. Saline

\* Wash For Ouchterlony Slides

For 2 liters:

200 ml 5% Sodium Bicarbonate  
1800 ml Normal Saline \*

1. Add the above <sup>together</sup> and it should  
make exactly 2 liters.
2. Now ready to use as wash  
for Ouchterlony slides.

\* Normal Saline = 9gms NaCl / 1 l. d. H<sub>2</sub>O.